

## Important Notice Regarding the Observation of Fixed Samples

Confocal microscope and STED microscope are used for both live and fixed specimen imagings. When you use these microscopes for fixed specimens, please be aware to use the following points.

1. Please seal the fixed cells and tissues in the slide glass tightly and be careful not to touch the objective lens with fixatives.
2. In case the fixatives touch the objective lens or microscope, please wash them immediately with water and alcohol. After that, please be sure to let us know it to the imaging core office (Ms. Asahara, extension: 28239, mail address: [asahara.masumi@mail.u-tokyo.ac.jp](mailto:asahara.masumi@mail.u-tokyo.ac.jp)).

Satoru Kondo  
IRCN Imaging Core